



## DSX Application Software Ver. 3.1.9

### Release Note

System Configuration	
OS	Microsoft Windows7 Professional 32bit/64bit SP1 Microsoft Windows10 Pro (Version 1511) 64bit Microsoft Windows10 Pro (Version 1607) 64bit Microsoft Windows10 Pro (Version 1703) 64bit Microsoft Windows10 Pro / Pro for Workstations (Version 1709) 64bit Microsoft Windows10 Pro / Pro for Workstations (Version 1803) 64bit Microsoft Windows10 Pro / Pro for Workstations (Version 1809) 64bit Japanese or English
.NET Framework	4.6.1 / 4.7 / 4.7.1 / 4.7.2
CPU	Intel Xeon®E3-1230 Quad-Core3.2GHz or more
RAM	4GB or more
HDD	500G or more ※The capacity to save images, etc. by the DSX application software depends on the free space in HDD.
Graphic	Equivalent to NVIDIA Quadro 600 1GB or more.
Optical Drive	DVD drive (The offline software can be installed to CD drive.)
Monitor Resolution	1920 x 1080



## 1. Changes from the previous version

### 1-1 Following features were supported.

No	Function	Description
1	To support newer Windows OS	Updated to support Windows10 Pro/ Pro for Workstations (Version1809) 64bit.

### 1-2 Following phenomena were corrected

No	Function	Description
1	Start-up	Fixed a bug that the application can not be started at the next time when the system time of the controller is changed before the application start time while the application is running.
2	Image List	Fixed a bug that the image list could not be closed when deleting an image opened from the image list from an external application such as Explorer.
3	Image List	Fixed a bug that the application is terminated when deleting a folder containing an opened image from an external application such as Explorer.
4	Image List	Fixed a bug that the application may be terminated when you select a large number of files displayed in the image list and press the delete button.
5	GUI related	In case of selecting Korean language, for the message displayed when the number of 3D shooting steps is 1000 or more, the number of steps shown as limit value 100 is corrected to 1000.



## 2. Compatibility table of OLYMPUS Stream

When used in combination with OLYMPUS Stream 2.1/2.2/2.3, please upgrade the DSX application software to 3.1.3 and after.

When used in combination with "OLYMPUS Stream 1.8 - 1.9.3" and "DSX application software before 3.1.2", after version up to the "DSX application software 3.1.3/3.1.3a", please version up to the "OLYMPUS Stream 1.9.4".

		OLYMPUS Stream					
		1.7	1.8~1.9.3	1.9.4	2.1	2.2	2.3
DSX-BSW	Before 3.1.2	✓	✓	✓	-	-	-
	3.1.3-3.1.4	✓	-	✓	✓	✓	✓
	3.1.5-3.1.9	✓	-	-	✓	✓	✓

## 3. Restrictions of Specifications

### 3-1 Restrictions remaining from the previous version (Ver. 3.1.8)

No	Function	Description	Countermeasures
1	Zoom Change	If the image acquired with the digital zoom magnification (the status when the [D] icon is displayed on the slider bar for specifying the magnification) is displayed in [vertical shrinking] or [horizontal shrinking] in [dual screen], the position to display the live image is shifted to the upper-left.	If comparing images in the dual screen, do not use the image acquired with the digital zoom magnification.
2	Z Movement	If Z movement is performed to the Z limit position while setting [Reference Z position] same as [Z limit position], only the message indicating that it reached [Reference Z Position] is displayed, but the message indicating it reached [Z Limit Position] is not displayed.	This is only a display issue and there is no functional problem.
3	Automatic Stitching	After completing the automatic stitching, the [Stop] button is not returned to the [Acquisition] button infrequently. If the [Stop] button is pressed during this status, the [Acquisition] button cannot be pressed.	This problem is solved by following procedures. (1) Close the automatic stitching screen. (2) Display the [Live Stitching] screen. (3) Press the [Acquisition] button.

4	HDR	If the HDR image is acquired while the live display stops, even though the fixed parameters for HDR detail settings is set to OFF, the automatic adjustment process of parameters is not executed.	Enable the live display before acquiring the image.
5	Display of Image Information	If the image acquired by using Ver1.1.5 and enabling HDR (Halation Removal) is opened with the application of this version, HDR=OFF is displayed in the image information display area.	This is only a display issue, and the opened image contains the information of HDR(Halation Removal) : ON.
6	GUI related	The top menu of the software (area where [Acquisition], [Measurement] or [Report] buttons are placed at the uppermost area of the screen) is not displayed in some cases.	Press the Windows key on the keyboard, and click the DSX button on the task bar displayed.
7		If [Acquisition], [Measurement] or [Report] screens are changed consecutively, it may take some time to change the screen.	Wait until the screen changes.
8		After the Z escape value is entered, the dotted line is displayed vertically.	This is only a display issue and there is no functional problem.
9		If the image in 3D display is opened while the multi-screen is displayed, the image is not displayed correctly in some cases.	Move the mouse cursor on the image display area, and then move it slightly.
10	GUI related	If [Acquisition] is selected from [Measurement] in Top Menu very quickly, the screen may not be displayed correctly.	Select [Acquisition] from [Measurement] in Top Menu again.
11		Immediately after the screen for registering users in Multi Preview is closed by clicking [Cancel], the acquisition screen may not be displayed correctly.	Click the [Measurement] tab or the [Acquisition] tab to display the screen correctly.
12		At re-login, when the message related to the anti-halation unit is displayed, the drawing of the background is not completed.	The message only appears during drawing the background. There is no functional problem.
13	Touch Panel	When moving the stage by instructing on the live image with the touch panel, if the next instruction is made before completing the last stage movement, the message indicating "The last operation may not have been executed." may appear.	After selecting [Image Correction]-[HDR], cancel the selection and operate the stage.

14	Observation setting	If the reproduction of the condition to lower the zoom magnification is repeated twice continuously, the control of the entire application will be Disabled.	Change the objective lens.
15	Step Measurement	If the image is opened in the order of 3D→2D→3D, the menu shows Plane Measurement, but the icon shows Step Measurement.	Nothing special
16	Report	If the size of the image opened in logo is large, it takes some time to save the report. ※ Several minutes for several MB.	Please wait for a while.
17		If the image is zoomed during saving the report, the internal error occurs.	Close the error message
18	Traveling acquisition	If the [Stop] button is clicked when the traveling acquisition of the first point is completed, the acquired image may not be displayed.	Go to the measurement screen. The acquired image is displayed there. Save it if necessary.
19	Video acquisition	After Video acquisition was finished in time out, Live observation and Video acquisition cannot be started.	Please restart application.